velleman®

CAMCOLMS

DIGITAL MICROSCOPE
DIGITALE MICROSCOOP
MICROSCOPE NUMÉRIQUE
MICROSCOPIO DIGITAL
DIGITALES MIKROSKOP

USER MANUAL
GEBRUIKERSHANDLEIDING
NOTICE D'EMPLOI
MANUAL DEL USUARIO
BEDIENUNGSANLEITUNG

1. Introduction

To all residents of the European Union

Important environmental information about this product

This symbol on the device or the package indicates that disposal of the device after its lifecycle could harm the environment. Do not dispose of the unit (or batteries) as unsorted municipal waste; it should be taken to a specialized company for recycling. This device should be returned to your distributor or to a local recycling service. Respect the local environmental rules.

If in doubt, contact your local waste disposal authorities.

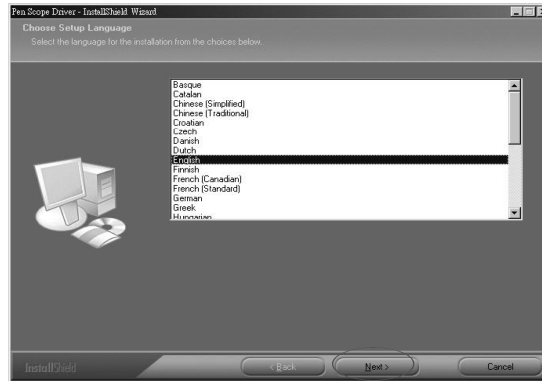
Thank you for choosing Velleman! Please read the manual thoroughly before bringing this device into service. If the device was damaged in transit, do not install or use it and contact your dealer.

2. System Requirements

- Pentium® MMX™ 200MHz with 32MB RAM or higher
- Windows® 98/ME/2000/XP/Vista
- free USB 2.0 or USB 1.1 port
- DirectX® 8.0 or higher

3. Installing the Driver

- Make sure the microscope is not connected to the computer. Insert the included CD-ROM into the CD-ROM drive of your computer. Select your language and click **[Next]**.

- Click **[Next]**. After the driver installation is completed, you will be asked to restart your computer. This is not necessary but recommended.

4. Installing the Microscope

- Connect the microscope to the computer through a USB port. A wizard will pop up. Choose not to search for software and click **[Next]**.

- Your computer automatically finds the connected microscope. Click **[Install the software automatically (Recommended)]** and click **[Next]** to continue the installation.

- Wait for the installation procedure to complete. Click **[Continue Anyway]** when following window pops up.

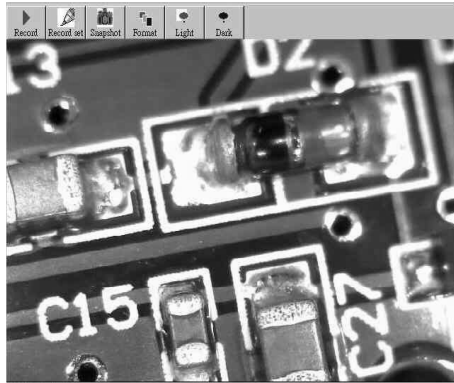
- Click **[Finish]** to complete your installation. Note that the microscope must only be connected to the USB port at time of installation. The computer will not recognize the microscope when connected to a different USB port.


5. Operating Instructions

- Place the microscope perpendicularly with and near the object to be viewed. Turn the adjustment wheel slowly to obtain as clear an image as possible.

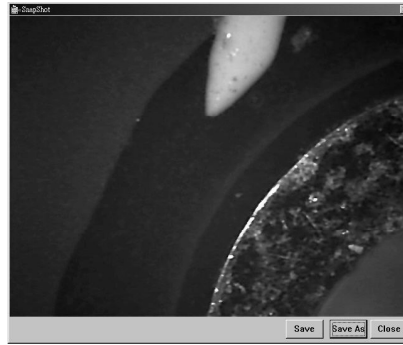
Click **[Snapshot]** to capture the screen. This screen capture can be saved as a JPEG or BMP file (JPG default). To do so, click **[Save]**. The file will be stored automatically in **[My Documents\PEN SCOPE\SNAPSHOT]** according to the following format: 070517001.JPG (07 = year, 05 = month, 17 = day, 001 = snapshot number, JPG = file format). Click **[Save As]** to configure the file name.

Click **[Record]** to make a recording.


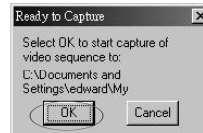
Click **[Record Set]** to make a recording according to your own specifications. Set the recording time and file size limit, click **[Start]** and **[OK]** to start recording. Click **[Save As]** to save the recording. The file will be stored automatically in **[My Documents\PEN SCOPE\RECORDING]** according to the following format: 070517001.AVI (07 = year, 05 = month, 17 = day, 001 = recording number, AVI = file format).

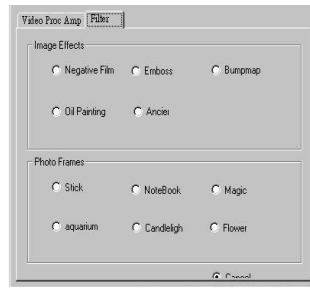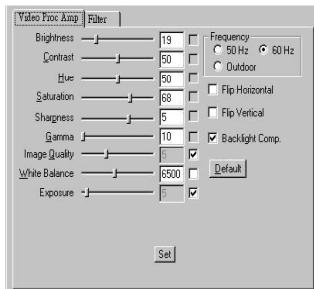Click **[Format]** to configure your screen settings or to add effects.

Click **[Light]** or **[Dark]** to switch the microscope on or off.

6. Technical Specifications

Image Sensor	CMOS
Number of Pixels	300,000
Effective Pixels	645 (H) x 485 (V)
Input Interface	USB 1.1 or 2.0
Magnification Rate	50x / 200x
Min. Illumination	10 ~ 20 lux / F2.8 – 5,600K
S/N Ratio	42dB
Cable Length	1.2m
White Balance	auto
White Balance Range	2,800 ~ 11,000K
Iris Aperture	AES
Shutter	1/30 ~ 1/15,750 sec
Picture Format	JPEG / BMP
Power Consumption	65 ~ 100mA
Dimensions	Ø 38 x 165mm
Weight	128g

Use this device with original accessories only. Velleman nv cannot be held responsible in the event of damage, injury or data loss resulted from (incorrect) use of this device.

For more info concerning this product, please visit our website www.velleman.eu.
The information in this manual is subject to change without prior notice.

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