velleman®

DVM8681

pH METER
pH-METER
pH-MÈTRE
pH-METRO
pH-MESSGERÄT

USER MANUAL
GEBRUIKERSHANDLEIDING
NOTICE D'EMPLOI
MANUAL DEL USUARIO
BEDIENUNGSANLEITUNG

1. Introduction

To all residents of the European Union

Important environmental information about this product

This symbol on the device or the package indicates that disposal of the device after its lifecycle could harm the environment.

Do not dispose of the unit (or batteries) as unsorted municipal waste; it should be taken to a specialized company for recycling.

This device should be returned to your distributor or to a local recycling service.


Respect the local environmental rules.

If in doubt, contact your local waste disposal authorities.

Thank you for choosing Velleman! Contents: 1 x pH meter, 4 x LR44 button cell and this manual. Please read the manual thoroughly before bringing this device into service. If the device was damaged in transit, don't install or use it and contact your dealer.



2. Battery Installation

- Loosen the two screws holding the battery cover. Do not discard the small washer.
- Insert 4 x LR44 button cell batteries respecting the polarity.
- Close the battery compartment.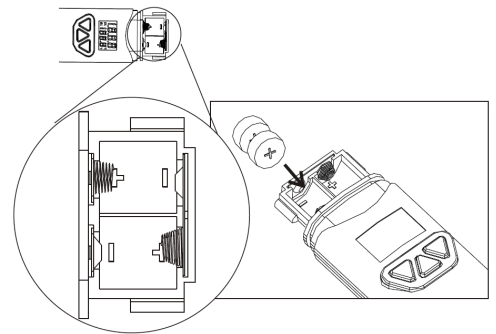

3. Description


1. measured pH reading
2. CAL – calibration mode
3. ATC – automatic temperature compensation
4. temperature reading
5. HLD – data hold
6. PWR – power button
7. CAL – calibration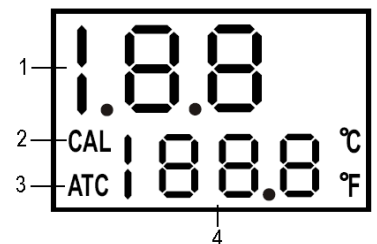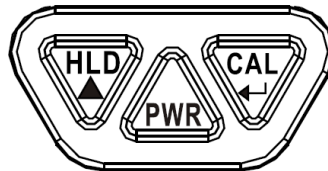


4. Operation

a. General Operation

- Remove the bottom and the inner cap to expose the electrode. It is normal to find some whitish crystals on the inner cap and electrode. **Keep the sponge in the cap wet to keep the electrode in a good storage condition.**
- Dip the electrode into the test solution. Press PWR and stir the pH meter to get a stable reading.
- A small dot will be flashing while the meter is in measuring mode. The temperature (in °C or °F) will also be displayed.
- Press HLD to freeze the current reading. Note that the dot will stop flashing. Press HLD again to release the reading.
- Turn off the meter by pressing PWR. Cover the electrode with the cap and store the meter at an ambient temperature between 0°C ~ 50°C (32°F ~ 122°F). Refer to “**d. Calibrating your Meter**” to perform the 3-points calibration.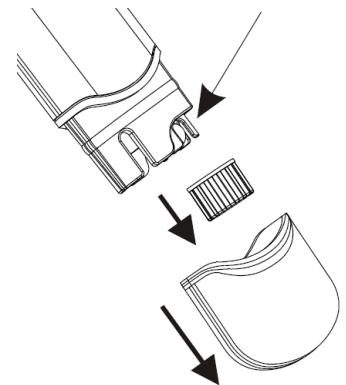


b. Auto Power-Off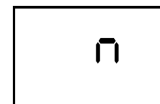

- This meter will switch off after approximately 20 minutes of inactivity.
- Disable this function by holding HLD and PWR simultaneously pressed before power-on. An “n” will briefly appear onto the display. The function has been disabled. **The function will be reactivated after every power-off.**


c. Automatic Temperature Compensation (ATC)

- Since pH values are temperature dependent, pH applications require some form of temperature compensation to ensure standardized pH values. This meter automatically corrects the pH values based on the temperature of the measured solution.
- Hold CAL and PWR simultaneously pressed before power-on. “C” or “F” will appear on the display. Select your temperature unit with HLD and press CAL to save.
- “SA” will briefly appear on the display to indicate the selection has been saved. **This selection will be saved after power-off.**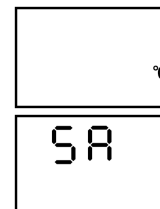


d. Calibrating your Meter

- Calibration is necessary and should be done regularly, even on a daily basis when using the meter often.
- Switch on your pH meter.
- Place the electrode into the pH 7 test solution first and into the pH 4 or pH 10 afterwards (test solution available at your local chemist’s).
- Press CAL to enter the calibration mode. “CA” will briefly appear on the display. After a few seconds “CAL” and the pH value will appear on the display indicating the meter has been calibrated.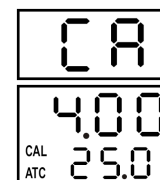

- The meter will display “En” if the calibration was not done correctly or if the probe could not detect the test solution in the voltage range below. In this case, recalibrate as per above.

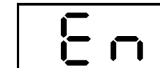| | |
|----------|-----------------|
| pH 4.00 | 97mV....250mV |
| pH 7.00 | -60mV....60mV |
| pH 10.00 | -250mV....-97mV |

- If the meter successfully recognizes the test solution, the solution pH value (4, 7 or 10) will appear on the display. If the displayed value differs slightly from the supposed value, e.g. 4.01 instead of 4.00, press HLD to change the value. The adjustable calibration range for pH 4 is 3.50 ~ 4.50, for pH 7 is 6.50 ~ 7.50 and for pH 10 is 9.50 ~ 10.50.
- A stable value will automatically be saved.

5. Cleaning and Maintenance

- Avoid touching the probe with bare hands and keep it wet by moistening the sponge in the cap.
- Rinse the probe with tap water or better, de-ionized water after each measurement.
- Remove the batteries from the meter if not used for a long time.

6. Troubleshooting

- **No display after power-on**
Check the batteries. Replace them if necessary (see “2. Battery Installation”).
- **Slow response**
Clean the probe thoroughly with tap water and rinse with de-ionized water afterwards.

- **Display indicates “- - -”**
The pH value is out of the pH range (pH 0 ~ pH 14). The solution is either too acid or too alkaline.
- **The displayed pH value is unstable**
The probe is exposed. Dip it in the solution to be measured.

7. Technical Specifications

| | |
|---------------------------|---|
| pH Range | pH 0 ~ pH 14 |
| Accuracy | pH 0.2 |
| Resolution | pH 0.1 |
| Temperature Accuracy | ± 1°C |
| Temperature Resolution | 0.5°C |
| Operating Temperature | 0°C ~ 50°C (32°F ~ 122°F) |
| Operating Humidity | 0 ~ 80% RH |
| ATC | yes |
| Self-Calibration | yes |
| Dimensions | 150 x 36 x 25mm |
| Calibration Test Solution | standard USA buffers, standard NIST buffers, standard DIN buffers |

Use this device with original accessories only. Velleman nv cannot be held responsible in the event of damage or injury resulted from (incorrect) use of this device.

The Velleman nv warranty covers normal operation and does not cover battery, misuse, abuse, alteration, tampering, neglect, improper maintenance or damage resulting from leaking batteries. Proof of purchase is required for warranty repairs. Warranty is void if the meter has been opened.

For warranty and service, please contact Velleman Components nv (www.velleman.eu).

For more info concerning this product, please visit our website www.velleman.eu.

The information in this manual is subject to change without prior notice.

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